

## Industry Application

## Speeding Up the Analysis of Dirty Water

Severn Trent Laboratories (STL) carry out analysis of dirty water samples including trade and sewage effluents, leachates, and contaminated ground water for organic parameters such as pesticides, phenols, and PAHs. One of STL's routine analytical methods involves the analysis of US EPA 8270 semi-volatile compounds (SVOCs) using Agilent GC-MSD systems operating in scan mode. In addition to providing results for about 100 target compounds, STL also produces a list of tentatively identified compounds (TICs) for each sample, together with the estimated concentrations. STL receives several thousand samples per year for this type of analysis.

Due to the complex nature of many of the chromatograms obtained in this process, the production of the TIC report is a highly skilled and time-consuming process. A software solution based around mass spectral deconvolution was therefore of interest.

STL needed a software solution that could bring a greater degree of automation to this process with deconvolution of samples, automatic search for compounds in the NIST database, and generation of a report with corresponding peak areas from which they could extract

semi-quantitative information. Intelligent Compound Recognition for MS Search powered by ACD/IntelliXtract (ACD/IXCR) performs all of these steps, and with reduced operator involvement.

After implementing ACD/IXCR and working with the software for a few months, STL found that deconvolution using the IntelliXtract algorithm was able to identify a wider range of compounds than by using the GC-MSD process alone. The software automatically searches the NIST database for similar mass spectral patterns (based on reconstructed pure component spectra) and produces a clear and concise report including names of identified compounds (with a match score) and the area under the peak. Importantly for STL, the report can be automatically exported into standard software where concentrations can be calculated.

Despite the software being very different to anything STL had previously used, they found it fairly intuitive to use, and the results were very comparable to other solutions they had evaluated. The real benefits for STL have come in the software's ability to help reduce analytical time, and therefore increase capacity for this type of analysis.



To learn more about how ACD/IXCR  
can work in your organization, visit  
[www.acdlabs.com/ixcr/](http://www.acdlabs.com/ixcr/) or email [info@acdlabs.com](mailto:info@acdlabs.com)

Visionary Software



Advancing Research